

QuickStart Setup Guide for SimplicityCam Microscope Adapter

REQUIRED COMPONENTS:

- 1) SimplicityCam Model: 5z, 5zAF or 8zAF.
- 2) Recordex Microscope Adapter (SC5zMA).
- 3) Standard microscope with 23mm eyepiece.
- 4) PC and USB Cable.

1. Make USB connection

Note, the SimplicityCam drivers will install when the camera is plugged in to the computer. No driver pre-installation is required. Simply attach the camera and wait a few minutes for Windows or OSX to install the drivers.



2. Remove the eyepiece from your microscope and replace with the SC5zMA adapter.

After replacing the eyepiece you can continue to use your microscope as normal. However, now it will also work with your SimplicityCam document camera.



3. Position the SimplicityCam directly over the microscope eyepiece.

Note, for detailed instructions view the training video located here:

<https://youtu.be/rRh15zl2ykM?list=PLJPmvEnrFMlixwwXoz4CSAW1QyREXX08K>



4. Go to www.recordexusa.com for more info

Visit our web site to download our doc cam software, XSight, to begin capturing images and video from your microscope using your SimplicityCam.

Copyright © 2016 RecordexUSA, Inc. All Rights Reserved.

Mac and Mac OSX are trademarks of Apple Inc., registered in the U.S. and other countries. Windows is a registered trademark of Microsoft Corporation in the United States and other countries. Recordex XSight software and Recordex SimplicityCam document cameras include a license to Pathway Innovations and Technologies, Inc.'s U.S. Patent No. 8,508,751.

recordex[™]